Considerations When Selecting Conical Tubes for Centrifugation Applications

Stephanie Carter, Robert Scott, Wenxiao Lu, Marci Moore, Joe Granchelli, Cindy Neeley

Key Words

Conical centrifuge tubes, centrifuge rotors and adaptors, RCF, RPM

Background

Laboratory centrifuges are common, everyday instruments. While centrifuge usage is widespread, the selection of the proper centrifuge and rotor system, along with the proper centrifuge tube consumable, can often be a challenging experience. There are several factors to consider when selecting the correct conical tube for centrifugation. These include the relative centrifugal force required, the fit of the particular tube in the rotor, sample volume, and the compatibility of the sample to the tube material. Understanding the requirements of a centrifugation application before selecting a conical tube can prevent sample leakage or loss, allow for easy sample recovery, and reduce the risk for potential damage to the centrifuge and rotor.



Factors to consider when choosing a conical centrifuge tube

A. RPM vs. RCF

All centrifuge tubes have a maximum speed rating determined by the manufacturer. Vessels used at speeds higher than the recommended rating can fail, resulting in sample loss and potential damage to the centrifuge and rotor. Most protocols specify speed in either revolutions per minute (RPM) or relative centrifugal force (RCF). It is crucial to understand the difference between RPM and RCF. The rotor revolves at the specified RPM and the force applied to the contents is dependent on the rotor's radius, with larger radii applying greater force on the sample. RCF represents the gravitational force being applied to the sample. It determines the centrifugation outcome independent of rotor size. RCF is measured in force x gravity, or g-force, and is more relevant to the actual impact and outcome of the centrifugation than RPM. Users should verify the required RCF by their specific applications to ensure that the required centrifugation force does not exceed the manufacturer's specified g-force rating.



B. Fit of the conical tube in the centrifuge rotor

It is important to note that the centrifuge tube performance largely depends on how well the tube fits in the rotor (or rotor adapter). For optimal tube performance it is imperative that there is contact between the tube and the rotor/adapter so that centrifugal force can be distributed to the rotor rather than the tube. Forces exerted directly on the conical tube may cause stress lines, bulging, or cracking. For example, when using a swingout bucket rotor in a centrifuge unit, the direction of the g force centers on the conical bottom of the tube locking the tube in the center position of the rotor. In some cases, this does not allow the lower portion of the tube to make contact with the rotor/adapter. The centrifugal forces on the side wall of the conical tube cannot be transferred to the rotor and may cause damage to the lower portion of the tube (Figure 1).

When using a fixed angle rotor in a centrifuge unit, the side of the tube makes contact with the rotor/adapter and most of the centrifugal force exerted on the tube can be transferred to the rotor (Figure 2). Consequently, the conical tube in a fixed angle rotor can be spun at much higher RCF in comparison to a swing-out bucket rotor.



Figure 1. Swing-out bucket rotor and tube fit illustration. Centrifugal forces are exerted on the conical and lower portion of the tube. The imperfect fit of the lower portion of the tube to the rotor leaves a gap which may cause tube damage.





Table 1 and Table 2 display the recommended maximum RCF for the Thermo Scientific Nunc 15 mL and 50 mL conical tubes based on tests of 9 models of rotors in 3

types of centrifuge units. In most cases, a rotor adapter is needed to ensure good fitting of the tube to the rotor.

Thermo Scientific Centrifuge Model	Thermo Scientific Rotor Model	Thermo Scientific Adapter Catalog #	Maximum RCF (x g)
Legend [™] 1XR	BIOShield [™] 720	75003678	3500
Legend 1XR	Fiberlite F15-6x100y	75003095	24652*
Legend 1XR	TX-400	75003682	4696*
Legend 1XR	TX-200	75003771	5580*
Multifuge [™] X3R	TX-750	75003639	4816*
Multifuge X3R	BIOShield 1000A	75003642	5000
Multifuge X3R	Fiberlite F15-8x50cy	010-0378-06	24446*
Multifuge X3R	BIOLiner	75003673	2739*
Evolution™	Fiberlite F13-14x50cy	010-0378-06	15000
LYNX 6000	Fiberlite A21-24x15c	not necessary	18500
LYNX 6000	BIOFlex HC	75007306	6500
LYNX 6000	BIOFlex HS	75005394	7000

Table 1. Recommended Maximum RCF for Thermo Scientific[™] Nunc[™] 15 mL Conical Tubes (Catalog # 339650, 339651)

* Maximum speed for the centrifuge/rotor system is reached.

Thermo Scientific Centrifuge Model	Thermo Scientific Rotor Model	Thermo Scientific Adapter Catalog #	Maximum RCF (x g)
Legend 1XR	BIOShield 720	75003677	6500
Legend 1XR	Fiberlite F15-6x100y	75003103	23500
Legend 1XR	TX-400	75003683	4696*
Legend 1XR	TX-200	75003803	5580*
Multifuge X3R	TX-750	75003638	4816*
Multifuge X3R	TX-750	75003824	4816*
Multifuge X3R	BIOShield 1000A	75003643	6500
Multifuge X3R	Fiberlite F15-8x50cy	not necessary	20000
Multifuge X3R	BIOLiner	75003674	2739*
Evolution	Fiberlite F13-14x50cy	not necessary	15000
LYNX 6000	Fiberlite F14-14x50cy	not necessary	13000
LYNX 6000	BIOFlex HC	75003674	7068*
LYNX 6000	BIOFlex HS	75005393	10025*

* Maximum speed for the centrifuge/rotor system is reached.

C. Chemical compatibility

In addition to the maximum RCF, when evaluating centrifuge tubes it is important to make sure that the sample components will not harm the plastic tube material. Chemicals can affect the strength, flexibility, surface texture, color, and shape of the plastic. Chemical resistance is influenced by temperature, duration and frequency of exposure, chemical concentration, and centrifugal force. Physical and chemical changes which may be caused by chemical exposure include:

- Absorption of solvents, resulting in softening or swelling of the plastic
- Stress-cracking of the plastic
- Permeation of solvent in the sample through the plastic
- Dissolution of polymer in the sample

The Nunc conical centrifuge tubes are made of premier quality polypropylene. Table 3 displays information on chemical compatibility of common reagents to polypropylene centrifuge tubes.

D. Sample volume

Before selecting a centrifuge tube size, the desired sample volume should be determined. The sample volume is limited to the available rotors in the laboratory and the protocol being used. Generally, a centrifuge tube should be filled at least 75%. In some cases, such as with ultracentrifuge tubes, it is required that tubes are completely filled to prevent failure. Using centrifuge tubes less than half filled to capacity can lead to high levels of material stress and can result in tube failure. If the available sample volume is smaller than the available rotor capacity, adapters are available for most rotors from the manufacturer (Table 1, Table 2). Adapters allow a smaller-volume tube to be used at the appropriate fill volume. This is essential when the recommended maximum RCF is going to be applied for centrifugation.

Classification	Chemical	Rating
Alcohols	Butanol, pure	S
	Ethanol, 100%	S
	Isopropanol, 100%	S
	Methanol, 15%	S
	Methanol, 98% x-750	S1
	Propanol, 100%	S
Cryopreservation Agents	Dimethyl sulfoxide (DMSO), pure	S
	Glycerol	S
Detergents	Sodium Dodecyl Sulfate (SDS), pure	S
	Triton X-100, pure	S
	Tween-20	S
Fixatives	Acetone, 50%	S
	Acetone, pure	Μ
	Formaldehyde, 10%	S1
	Formaldehyde, 30%	Μ
	Formalin, 10%	S1
	Formalin, 30%	Μ
	Gluteraldehyde, pure	S
	Paraformaldehyde, pure	Μ
Other	Beta- Mercaptoethanol, pure	Μ
	Cell culture media and sera	S
	EDTA, pure	S
	RNAzol	S
	Trypsin	S
	Toluene, pure	U
	Xylene, pure	U

Key:

S = Satisfactory

S1 = Satisfactory, may cause discoloration.

M = Marginal; may be satisfactory for use in a centrifuge, depending on length of exposure and speed. Testing under operating conditions is suggested before actual run.

U = Unsatisfactory; not recommended.

Summary

As in most endeavors, gathering the correct information ahead of time is essential. In this study, we demonstrated the important factors to consider when selecting the conical tubes for centrifugation so that sample leakage or loss is prevented. Choosing the correct centrifuge consumables also ensures proper execution of centrifugation, and reduces the risk for potential damage to the centrifuge and rotor. The factors we discussed in this article include the RCF required by the protocol, the fit of the particular conical tube in the rotor, sample volume, and the chemical compatibility of the sample to the tube material. Users should verify the maximum RCF of the centrifuge tube to ensure that the required speed does not exceed the manufacturer's rating. Here, we recommended the maximum RCF of the Nunc 15 mL and 50 mL conical tubes in several Thermo Scientific centrifuge and rotor systems while taking these factors into consideration.

References

Goodman, Tammy (Thermo Scientific). Selecting Centrifuge Consumables. American Laboratory April 2009.

Thermo Scientific, NALGENE Centrifuge Tubes and Bottles Your Complete Guide to Centrifuge Ware, 2001

Ordering Information

Catalog Number	339650	339651	339652	339653
Volume, mL	15	15	50	50
Packaging	Bulk	Racked	Bulk	Racked
Material	PP	PP	PP	PP
Units per pack/case	50/500	25/500	25/500	25/300

All centrifuge tubes are Sterile - SAL 10⁻⁶, Leak proof, USP Class VI, Non-pyrogenic, Non-cytotoxic and RNase/DNase Free.

thermoscientific.com

© 2014 Thermo Fisher Scientific Inc. All rights reserved. All (other) trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries. Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details.

ANZ: Australia: 1300 735 292, New Zealand: 0800 933 966; Asia: China Toll-free: 800-810-5118 or 400-650-5118; India: +91 22 6716 2200, India Toll-free: 1 800 22 8374; Japan: +81-3-5826-1616; Other Asian countries: 65 68729717 Europe: Austria: +43 1 801 40 0; Belgium: +32 2 482 30 30; Denmark: +45 4631 2000; France: +33 2 2803 2180; Germany: +49 6184 90 6000, Germany Toll-free: 0800 1-536 376; Italy: +39 02 95059 554; Netherlands: +31 76 571 4440; Nordic/Baltic countries: +358 9 329 10200; Russia/CIS: +7 (812) 703 42 15; Spain/Portugal: +34 93 223 09 18; Switzerland: +41 4454 12 22; UK/Ireland: +44 870 609 9203 North America: USA/Canada +1 585 586 8800; USA Toll-free: 800 625 4327 South America: USA sales support: +1 585 899 7198 Countries not listed: +49 6184 90 6000 or +33 2 2803 2000

Thermo s c I E N T I F I C A Thermo Fisher Scientific Brand