MP Biomedicals

Case study - Lactic Acid Bacteria (LAB)

Comparaison of three methods for determination of protein concentration in lactic acid bacteria for proteomics studies.

Ibrahim Mehmeti, Fadime Kiran, and Ozlem Osmanagaoglu Afr. J. Biotechnol. 2011. Vol 10.

Overview

- Keywords: FastPrep®, Sonication, centrifugation, lactic acid bacteria (LAB)
- Aim of the study: Development of an optimized protein extraction protocol
- Application: One dimensional (1D) SDS-PAGE. Two dimensional (2D)-PAGE
- Sample type: Bacteria (E. faecalis, P. pentosaceus and L. lactis)
- Material: Sonicator, centrifuge, FastPrep® 120 instrument
- Buffers: 8M Urea, 2M Thiourea, 0,5% CHAPS, 10 mM DTT and 0,1% immobilized pH gradient (IPG)

Protocol and Parameters

- 1. Pellets were resuspended in 400 µL rehydration buffer containing 9M Urea, 2M Thiourea, 0,5% CHAPS, 10 mM DTT and 0,1% immobilized pH gradient (IPG).
- 2. Cells were lysed with acid washed glass beads (diameter of 212 to 300 μ m) using a FastPrep® 120 at speed 6 m/s for 3 x 45 sec. at 4°C. After each cycle the solution was kept on ice for 1 min.
- 3. Cell debris were removed by centrifugation at 14,000 x g for 10 min at 4°C.
- 4. Before analysis, the supernatant was kept at -20 °C.

Results

Six fold greater amount of protein was obtained with FastPrep® bead beater:

Method	Sonication	Centrifugation	FastPrep
E. faecalis V583	1.33 ± 0.01	1.25 ± 0.01	4.85 ± 0.05
L. lactis NIZO 0900	1.25 ± 0.02	1.32 ± 0.01	6.23 ± 0.06
P. pentosaceus OZF	1.32 ± 0.01	$1,16 \pm 0.02$	5.66 ± 0.04


Mean \pm SD of protein concentrations (μ g/ μ l) of each strain obtained by three different methods. Values are mean \pm S.D.(standard deviation) of results of three experiments.



www.mpbio.com


Results

Higher efficiency and higher quality in extracting proteins with FastPrep® method:

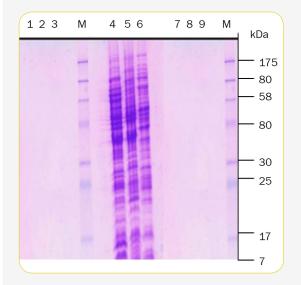

Representative Coomassie brilliant blue stained SDSPAGE illustrating the intracellular proteins of three representative strains of LAB. The lanes (1 to 9) contains extracts of P.pentosaceus OZF (lanes 1, 4, 7); E. faecalis V583 (lanes 2, 5, 8) and L. lactis NIZO 0900 (lanes 3, 6, 9) obtained by centrifugation (lanes 1 to 3), FastPrep® (lanes 4 to 6) and sonication (lanes 7 to 9); M, prestained broad range protein marker (Bio Labs).

FastPrep® method showed higher protein resolution and spot intensity of all proteins:

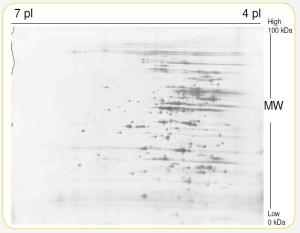

Representative 2D-PAGE illustrating the intracellular proteins of E. faecalis V583 at pl 4 to 7 after lysing cells by FastPrep®. See higher protein resolution and spot intensity.

Conclusion

- The results show that higher amount of proteins were obtained when the cells were lysed with a FastPrep®. **Six times higher protein concentration**, which is important for proteomics, was obtained with extraction by FastPrep®.
- SDS-PAGE gel images show that higher amounts of proteins are obtained only when proteins extracted by FastPrep® method.
- More than 400 protein, spots, with isoelectric points (pl) ranging from 4.0 to 7.0 and molecular weights (MW) from 0 to 100 kDa, were observed with 2D-PAGE analysis. In addition, proteins with less abundance and high molecular weights were resolved clearly and detected strongly on 2D gel when FastPrep® method was used.
- FastPrep® extraction method was an efficient and reliable method for lysing and/or extracting proteins of *LAB* for proteomic approach and reproducible amounts of bacterial proteins can successfully be extracted.
- Pictures were excellent enough to be used in alignment for statistical analyses and spots well-resolved for MALDI TOF analyses.

Successful sample preparation using the MP Biomedicals FastPrep® product line has been highlighted in thousands of scientific articles. To access articles and other materials, visit www.mpbio.com/FastPrepLibrary.



www.mpbio.com