

Optimized technology and ergonomic design!



macro Pipette Controller

FIRST CLASS · BRAND

The macro pipette controller from BRAND – comfortable and simple pipetting!

- Draws liquids up rapidly
- Easy, sensitive meniscus adjustment with the spring-loaded control lever
- Designed without corners and edges for convenient pipetting, even for long serial pipetting
- Four colors to choose from: Neutral gray, bold magenta, fresh green, or friendly blue, for better identification of different applications



At the www.brand.de site, you will find a large selection of volumetric and measuring pipettes, now also compliant with USP requirements.



Convenient operation without effort

The unique valve system allows for easy compression of the bellows. 50 ml of fluid can be drawn in within just 11 seconds. The spring loaded lever enables sensitive filling and delivery of liquids. The meniscus is easily adjustable.

Ergonomic design

The optimized design, the practical arrangement of functions, and the low weight of 125 g offer reliable operation during serial pipetting tasks, even for inexperienced users.



Broad area of applications

One single macro pipette controller covers the entire range of bulb and graduated pipettes from 0.1 to 200 ml. The conical silicone adapter offers a secure fit for the different diameters. The unit is fully autoclavable at 121 °C (2 bar) according to DIN EN 285. A hydrophobic membrane filter protects the system from liquid penetration.

Perk up the lab with color

Four different colors: gray, magenta, green, and blue.

Effortless Pipetting



Filling

Squeeze the suction bellows. It allows aspiration up to 50 ml with one compression. The fill rate is continuously variable with the pipetting lever.

Ordering Data

macro Pipette Controller, incl. spare membrane filter, 3 µm.
Pack of 1.

| Color | Cat. No. |
|---------|----------|
| gray | 262 00 |
| magenta | 262 03 |
| green | 262 01 |
| blue | 262 02 |

Spare parts

| Description | Pack of | Cat. No. |
|---|---------|----------|
| Membrane filter 3 µm (PP, PTFE), non-sterile | 1 | 260 52 |
| Membrane filter 3 µm (PP, PTFE), non-sterile | 10 | 260 56 |
| Adapter (Silicone), length 44 mm | 1 | 261 46 |
| Adapter support (PP), length 49 mm | 1 | 262 20 |
| Suction bellows (Silicone) with screw ring (PP) | 1 | 262 25 |
| Valve system (PP, PTFE, Silicone) | 1 | 261 28 |

BLAUBRAND® Pipetting Package



Items supplied:

- **1 macro pipette controller**, gray
- **6 BLAUBRAND® graduated pipettes, type 2**
3 pipettes 5 ml and 10 ml each, class AS, DE-M marking, with batch certificate
- **Useful product information**
BLAUBRAND® volumetric instruments, handling of pipettes
- **Handy plastic container**
Ideal for storage of pipettes up to 360 mm length

Cat. No. 260 07



Delivery

To discharge the liquid, simply press down the lever. The valve system is designed so that flow rate does not affect accuracy.



Blow-out function

When using blow-out pipettes, press the convenient blow-out bulb once the meniscus has come to rest in the jet. Pipetting has never been easier.

BRAND® and BLAUBRAND® are trademarks of BRAND GMBH + CO KG, Germany.

Our technical literature is intended to inform and advise our customers. However, the validity of general empirical values, and of results obtained under test conditions, for specific applications depends on many factors beyond our control. Please appreciate, therefore, that no claims can be derived from our advice. The user is responsible for checking the appropriateness of the product for any particular application.

Subject to technical modification without notice. Errors excepted.

BRAND GMBH + CO KG · P.O. Box 11 55 · 97861 Wertheim · Germany
Tel.: +49 9342 808-0 · Fax: +49 9342 808-98000 · E-Mail: info@brand.de · Internet: www.brand.de

